



ken-a-vision[®]
KNOWLEDGE THROUGH VISION

Microscope Replacement Heads

Instruction Manual

SC12MH

SC12DH

SC12BH

SC12CMH

SC12CBH



Figure 1

Place your 5/64 (1.98MM) Allen wrench in the Allen set screw located on the right side of the microscope. Figure (1)



Figure 2

Loosen screw carefully remove the existing head. Figure (2)



Figure 3

Locate your new replacement head. Be sure to remove plastic cap located on the base of the head. Figure (3)



Figure 4

Carefully place the replacement head in the cavity, align and tighten allen set screw. Figure (4)



ken-a-vision®
KNOWLEDGE THROUGH VISION

www.ken-a-vision.com
INS-SHEAD